



Instruments

Pitt County Sheriff's Office Forensics Services Unit
Issued by the Technical Leader

Effective
Date:
2018/04/01

Ver:
2

Drug Chemistry Technical Procedure for Polarized Light Microscopy

Page #:
1 of 9

1.0 Purpose - This procedure specifies the required elements for the preparation and use of microcrystalline reagents, and viewing suspected marijuana/ hashish/THC samples with the polarized microscope.

2.0 Scope - This procedure applies to all polarized light microscopy techniques used in the Drug Chemistry section of the Pitt County Sheriff's Office Forensic Services Unit.

3.0 Definitions

- **Quality control (QC) check** - Periodic confirmation of the reliability of equipment, instrumentation, and/or reagents.
- **Reference Material** – Material sufficiently homogeneous and stable with reference to specified properties, which has been established to be fit for its intended use in measurement or in examination of nominal properties.

4.0 Equipment, Materials and Reagents

4.1 Equipment

- *Leica* DM750P Polarized light microscope

4.2 Materials and Reagents

- Fume hood
- Eye protection
- Laboratory coat
- Gloves
- Balance
- Beakers or other glass vessels
- Graduated cylinder
- Glass stirring rod
- Reagent bottle(s)
- Microscope slides
- Objective centering screws
- Spatula
- Weigh boats or other weigh vessels
- Reference materials

5.0 Procedure

5.1 Standards and Controls - Quality control checks of all reagents shall consist of a negative check and a positive check. Both checks shall be acceptable according to the procedure listed for each reagent, and shall be recorded together as a quality control check in the Reagent Log.

5.1.1 Negative quality control checks shall be performed according to the procedure listed with no sample present.

All copies of this document are uncontrolled when printed.



Instruments
Pitt County Sheriff's Office Forensics Services Unit
Issued by the Technical Leader

Effective
Date:
2018/04/01

Ver:
2

Drug Chemistry Technical Procedure for Polarized Light Microscopy

Page #:
2 of 9

5.1.1.1 Acceptable results are: no crystal formation or no hairs present.

5.1.1.2 If crystals do form or if hairs are present, steps will be taken until no crystals are formed or no hairs are present. This includes retesting with a new microscope slide, re-cleaning any utensils used, or making a new reagent/obtaining new solvent.

5.1.2 Positive and negative quality control checks shall be performed at six month intervals according to the procedure listed for each reagent using the reference material listed. See each procedure for acceptable results.

5.1.2.1 The result of the quality control check shall be recorded in the Reagent Log according to the [Drug Chemistry Technical Procedure for Quality Assurance](#).

5.1.3 Microcrystalline reagents may be prepared in any amount provided that the component ratios are kept constant.

5.1.4 Microcrystalline reagents shall expire three years from date of preparation.

5.1.5 Reagent bottles shall be labeled and checked as described in the [Drug Chemistry Technical Procedure for Quality Assurance](#).

5.2 Operation of the Polarized Light Microscope

5.2.1 Switch on the light source. (Refer to the operator manual for location/description of specific parts mentioned below.)

5.2.2 Place the specimen slide on the stage.

5.2.3 Adjust the desired light intensity with the control lever.

5.2.4 Make sure the field diaphragm is open to the edge of the field view.

5.2.5 Focus with the coarse and fine adjustments for the desired objective.

5.2.6 Move the microscope slide around to view the entire specimen, adjusting the focus accordingly.

5.2.7 Push the filter in to view the specimen with polars crossed, or pull it out to view with uncrossed polars.

5.2.8 If the objective is changed:

5.2.8.1 Adjust the fine focus adjustment.



Instruments
Pitt County Sheriff's Office Forensics Services Unit
Issued by the Technical Leader

Effective
Date:
2018/04/01

Ver:
2

Drug Chemistry Technical Procedure for Polarized Light Microscopy

Page #:
3 of 9

5.2.8.2 Set the field diaphragm to just inside the field of view.

5.2.8.3 Adjust the aperture diaphragm for optimum contrast and resolution.

5.3 Application of Procedures on Evidence

5.3.1 Dry Sample Method

5.3.1.1 A small portion of sample shall be placed on a microscope slide and a drop of the reagent or solvent shall be mixed with the sample.

- Optional step indicated in some procedures uses an acid solution (B) which may be used to dissolve the sample before addition of the reagent solution (A).

5.3.1.2 Any crystals formed or hairs present shall be observed under non-filtered and/or filtered light.

5.3.2 Criteria for Establishing Matches

5.3.2.1 Test samples that have evaporated to dryness may not be used for evaluation of crystals.

5.3.2.2 Crystals are identified by morphology (i.e., shape).

5.3.2.3 Crystals or cystolithic hairs shall be compared to those of standards tested in the same manner unless frequently observed.

5.3.2.4 Final comparison shall be with actual crystals given by the standards listed.

5.4 Formulas for Preparing Reagents

5.4.1 5 % Mercuric Chloride

5.4.1.1 This reagent is used for heroin and caffeine.

5.4.1.2 5 % Mercuric Chloride (A)

5.4.1.2.1 Add 1.5 grams of mercuric chloride to 30 milliliters of distilled water (5% weight volume solution).

5.4.1.2.2 Suggested Lot number format:
year/month/day/MerCur5/Initials of preparer.

5.4.1.3 0.05 N Hydrochloric Acid (B)



5.4.1.3.1 Add 1 milliliter of concentrated hydrochloric acid to 250 milliliters of distilled water.

5.4.1.4 Application of Procedure on Evidence

5.4.1.4.1 Mix directly with a drop of 5% mercuric chloride (A).

5.4.1.4.2 Optional step: Dilute a small portion of the sample on a microscope slide in a drop of 0.05N hydrochloric acid (B) before mixing with a drop of the 5% mercuric chloride (A).

5.4.1.5 QC check: Heroin forms fans/dendrites

5.4.1.6 Results: Heroin forms fans/dendrites
Caffeine – dendrites, but longer and less dense than heroin dendrites.

5.4.2 Gold Chloride in 20 % Acetic Acid With Optional 0.05 N Hydrochloric Acid

5.4.2.1 This reagent is used for cocaine and phencyclidine.

5.4.2.2 Gold Chloride/20 % Acetic Acid (A)

5.4.2.2.1 Add 10 milliliters glacial acetic acid to 40 milliliters of water.

5.4.2.2.2 Dissolve 1.0 gram of gold chloride in the 50 milliliters of 20 % acetic acid, with stirring.

5.4.2.2.3 Suggested Lot number format:
year/month/day/GdCl₂0/Initials of preparer.

5.4.2.3 0.05 N Hydrochloric Acid (B)

5.4.2.3.1 Add 1 milliliter of concentrated hydrochloric acid to 250 milliliters of distilled water.

5.4.2.4 Application of Procedure on Evidence

5.4.2.4.1 Mix directly with a drop of gold chloride/20% acetic acid (A).

5.4.2.4.2 Optional step: Dilute a small portion of the sample on a microscope slide in a drop of 0.05N hydrochloric acid (B) before mixing with a drop of gold chloride/20% acetic acid (A).

5.4.2.5 QC Check: Cocaine forms cross shaped crystals.



Instruments
Pitt County Sheriff's Office Forensics Services Unit
Issued by the Technical Leader

**Effective
Date:**
2018/04/01

Ver:
2

Drug Chemistry Technical Procedure for Polarized Light Microscopy

Page #:
5 of 9

5.4.2.6 Results: Cocaine forms cross-shaped crystals.
Phencyclidine forms squares with diagonal markings, often elongated along one axis.

5.4.2.7 Limitations: This technique does not distinguish between the salt and free base forms or the d-/l- forms of cocaine. It can however distinguish cocaine from its diastereoisomers.

5.4.3 50 % Acetic Acid and Gold Chloride in 50 % Acetic Acid

5.4.3.1 This reagent is used for enantiomer determination for propoxyphene.

5.4.3.2 Gold Chloride in 50 % Acetic Acid (A)

5.4.3.2.1 Carefully add 20 milliliters glacial acetic acid to 20 milliliters of distilled water.

5.4.3.2.2 Mix the contents of a one gram ampoule of gold chloride into the 50 % acetic acid solution.

5.4.3.2.3 Suggested Lot number format:
year/month/day/GdCl150/Initials of preparer.

5.4.3.3 50 % Acetic Acid (B) – Carefully add 20 milliliters glacial acetic acid to 20 milliliters of distilled water.

5.4.3.4 Application of Procedure on Evidence - Mixed Crystal Testing (for enantiomer determination)

5.4.3.4.1 Mix a portion of the unknown sample directly with reagent (A), or dilute the sample on a microscope slide in a drop of 50 % acetic acid (B) before mixing with a drop of the Gold Chloride in 50 % acetic acid (A). Crystals formed are compared with crystals formed from the mixture(s) in the next step.

5.4.3.4.2 Mix a second portion of the unknown sample with an equal portion of the d- or l- isomer of propoxyphene. Mix directly with reagent (A), or dilute the sample on a microscope slide in a drop of 50 % acetic acid (B) before mixing with a drop of the Gold Chloride in 50 % acetic acid (A). Crystals formed are compared with crystals formed from the sample in the previous step.

5.4.3.4.3 Single isomer and mixed isomer crystals have different appearances. Compare the crystals formed from the straight unknown to the d- and l-mixed samples. Isomer determination



may be made when racemic crystals are formed by one of the known isomer additions.

5.4.3.4.4 QC Check: d,l-Propoxyphene forms small, curved, irregular needles almost at once.

5.4.3.4.5 Results: d,l-Propoxyphene forms small, curved, irregular needles almost at once.
Single isomer propoxyphene (d- and l-) give large, straight needles which are very slow to form.

5.4.4 Microscopic Examination of Marijuana/Hashish Using Chloroform

5.4.4.1 This reagent is used to identify plant particles of marijuana.

5.4.4.2 Chloroform shall be used.

5.4.4.3 Application of Procedure on Evidence

5.4.4.3.1 Place a small sample of suspected hashish or marijuana on a microscope slide and add a drop of chloroform to the material.

5.4.4.3.2 Observe under a relatively low magnification (approximately 10x).

5.4.4.3.3 QC Check: Marijuana or hashish cystolithic hairs look like bear claws.

5.4.4.3.4 Results: Marijuana or hashish cystolithic hairs look like bear claws.

5.5 New Microscopes

5.5.1 New microscopes shall be installed by an approved vendor according to the manufacturer's instructions, and documented in the Document Management (DM).

5.6 Maintenance

5.6.1 Polarizing microscopes shall be serviced yearly by an approved vendor, and service visits documented in DM.

5.6.2 When a microscope has been placed out of service (e.g., maintenance, malfunction, leaving the direct control of the Laboratory), correct operation shall be verified by a successful cocaine crystal QC check.



Instruments

Pitt County Sheriff's Office Forensics Services Unit
Issued by the Technical Leader

Effective
Date:
2018/04/01

Ver:
2

Drug Chemistry Technical Procedure for Polarized Light Microscopy

Page #:
7 of 9

5.6.3 Laboratory personnel shall examine the effect(s) if any, of a malfunction on analysis and implement the [Laboratory Procedure for Corrective Action](#) as required.

5.6.4 If the amount of light passing through the optics decreases significantly so that a sample cannot be seen, steps shall be taken to correct this. This may include, but is not limited to, checking the Kohler illumination parameters.

6.0 Limitations

6.1 Diluents may interfere with crystal formation in some cases. Extraction or solvent washes may be needed to remove unwanted components before microcrystalline reagents are used.

6.2 Concentration of samples may need to be increased or decreased to aid in crystal formation.

7.0 References

Leica DM750 User Manual (Downloaded .pdf from internet after purchase.)

Allen, A.C., Cooper, D.A., Kiser, W.O., Cottrell, R.C. "The Cocaine Diastereoisomers", *Journal of Forensic Sciences*, JFSCA, Vol. 26, No. 1, Jan. 1981, pp. 12-26.

Brinsko, K.M, Golemis, D., et. Al. *A modern compendium of microcrystal tests for illicit drugs and diverted pharmaceuticals*, McCrone Research Institute, Chicago; 155 pages; 2016.

"Particle Characterization by PLM. Part I No. Polarities." *Microscope*, Volume 30, Issue 3 (1982): 185-196.

Butler, William P. *Methods of Analysis for Alkaloids, Opiates, Marihuana, Barbiturates, and Miscellaneous Drug.*, Publication #341. Washington, D.C.: U.S. Treasury Department, Internal Revenue Service December, 1966: 55, 114.

Clarke, E.G.C., and R.G. Todd, eds. *Isolation and Identification of Drugs*. 1st Edition. London: Pharmaceutical Press, 1969: 135-141,801.

Fulton, C.C. *Modern Microcrystal Tests for Drugs – The Identification of Organic Compounds by Microcrystalline Chemistry*. New York: Wiley-Interscience, 1969.


Sobol, S.P. and R.A. Moore. *Analytical Manual*, J.W. Gunn, ed. Bureau of Narcotics and Dangerous Drugs, Washington, D.C.: U.S. Government Printing Office 0-506-836, 1970: 105.

Bureau of Narcotics and Dangerous Drugs Seminar, 1970.

Smith, F.P., ed. *Handbook of Forensic Drug Analysis*. Boston, Massachusetts: Elsevier Academic Press, 2005: 238.

ASTM International Standard E1968, 2011, "Standard Guide for Microcrystal Testing in Forensic Analysis of Cocaine." ASTM International: West Conshohocken, PA, 2011, www.astm.org.

All copies of this document are uncontrolled when printed.

	<i>Instruments</i> Pitt County Sheriff's Office Forensics Services Unit <i>Issued by the Technical Leader</i>	Effective Date: 2018/04/01	Ver: 2
Drug Chemistry Technical Procedure for Polarized Light Microscopy			Page #: 8 of 9

ASTM International Standard E-2125, 2007, "Standard Guide for Microcrystal Testing in the Forensic Analysis of Phencyclidine and Its Analogues." ASTM International: West Conshohocken, PA, 2007, www.astm.org.

8.0 Records

- Reagent Log(s)
- Microscope maintenance records in DM.



Instruments
Pitt County Sheriff's Office Forensics Services Unit
Issued by the Technical Leader

Effective
Date:
2018/04/01

Ver:
2

Drug Chemistry Technical Procedure for Polarized Light Microscopy

Page #:
9 of 9

REVISION HISTORY

CURRENT VERSION	EFFECTIVE DATE	SUMMARY OF CHANGES
1	2017/11/14	Original Document.
2	2018/04/01	Header – Added “Drug Chemistry” Scope – Updated “Illicit Drugs discipline” to “Drug Chemistry section”